

BCA Protein Assay Kit

Product code: K001

Kit Contents:

Kit Contents:		Storage			
Extraction Times		100	500	2500	
I	Reagent A	20ml	100 ml	500 ml	RT
II	Reagent B	1ml	5 ml	25 ml	RT
III	BSA(5 mg/ml)	0.4ml	2 ml	10 ml	-20°C

Storage:

The contents of the kit should be stored at specified conditions and warmed to room temperature before used for the best testing sensitivity.

Introduction

The BCA protein assay is a protein determination formulation based on bicinchoninic acid (BCA) for the colorimetric detection. This method combines the reduction of Cu_2^+ to Cu^+ by protein in an alkaline medium (the biuret reaction) and the soluble purple-colored reaction product from the complexing of Cu^+ and BCA. This purple-colored complex exhibits a maximum absorbance at 562 nm that is nearly linear with increasing protein concentrations over a broad working range (0-2 mg/ml). According to highly sensitivity and simple to use, the BCA protein assay is adopted by laboratories and companies, and becomes one of the prior protein quantitative methods as well as the Bradford assay.

The Fine BCA Protein Assay Kit has the characteristics of high sensitivity and light background, and the measured range can be up to 3 mg/ml.

Procedure

1. 96-Microplate Procedure

(1) Prepare the Working Reagent. Working reagent, clear liquid in the color of yellow green, is mixed with the

1

Wuhan Fine Biotech Co., Ltd.

B9 Bld, High-Tech Medical Devices Park, No. 818 Gaoxin Ave. East Lake High-Tech Development Zone, Wuhan, Hubei, China(430206)

Tel: (0086)027-87384275

Fax: (0086)027-87800889

www.fn-test.com

Reagent A and Reagent B at ratio of 50:1.

(2) Prepare the Standard Reagent. Dilute the 5 mg/ml BSA Standard into lateral concentrations of 0, 0.25, 0.5, 0.75, 1, 1.5, 2 mg/ml. The BSA dilutions can be frozen and stored at -20°C, and thawed and warmed to room temperature when used.

(3) Mark the wells with Standard or Test. Add at least 20 µl of the diluted Standard Reagent to each Standard well and add equal volume of testing samples each Test well.

(4) Add all the wells with 200 µl Working Reagent and incubate for 30 min at 37°C.

(5) Measure the absorbance at 562 nm on a plate reader. Subtract the average absorbance reading of at least 3 blank wells. Draw a standard curve by plotting absorbance reading for each BSA standard versus its concentration. Use the standard curve to determine the protein concentration of each unknown sample.

2. Test Tube Procedure

(1) Prepare the Working Reagent. Working reagent, clear liquid in the color of yellow green, is mixed with the Reagent A and Reagent B at ratio of 50:1.

(2) Prepare the Standard Reagent. Dilute the 5 mg/ml BSA Standard into lateral concentrations of 0, 0.25, 0.5, 0.75, 1, 1.5, 2 mg/ml. The BSA dilutions can be frozen and stored at -20°C, and thawed and warmed to room temperature when used.

(3) Mark the tubes with Standard or Test. Add at least 50 µl of the diluted Standard Reagent to each Standard tube and add equal volume of testing samples into each Test tube.

(4) Add all the tubes with 1 ml Working Reagent and incubate for 30 min at 37°C.

(5) Measure the absorbance at 562 nm with a spectrophotometer. Subtract the absorbance average reading of at least 3 blank tubes. Draw a standard curve by plotting absorbance reading for each BSA standard versus its concentration. Use the standard curve to determine the protein concentration of each unknown sample.


Wuhan Fine Biotech Co., Ltd.

B9 Bld, High-Tech Medical Devices Park, No. 818 Gaoxin Ave. East Lake High-Tech Development Zone, Wuhan, Hubei, China(430206)

Tel: (0086)027-87384275

Fax: (0086)027-87800889

www.fn-test.com

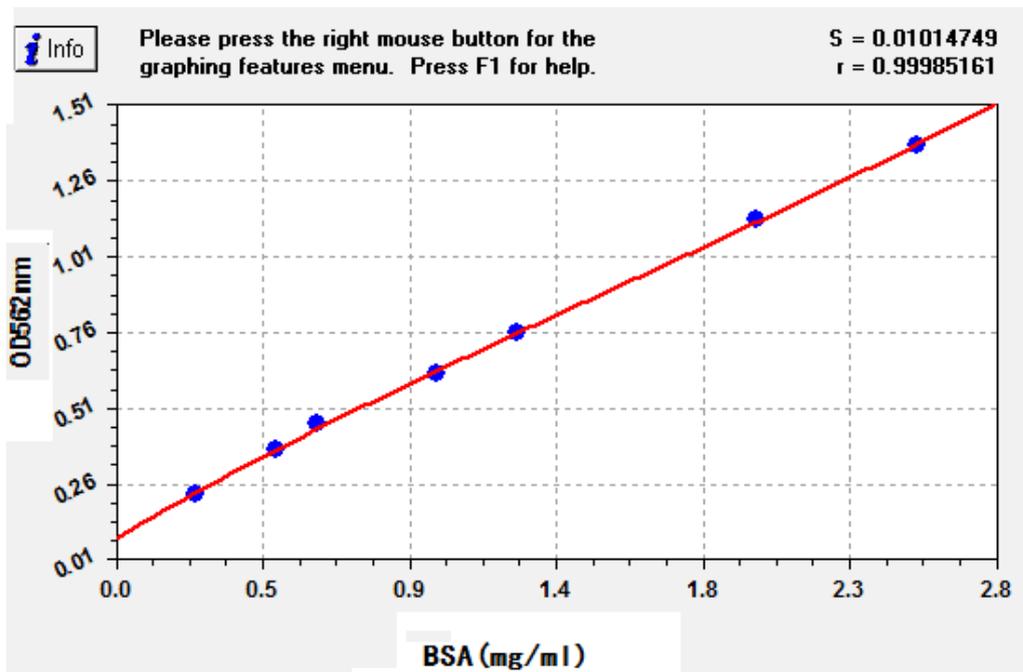

Additional Testing Information

- 1.If lacking time for incubation, the mixed reagents can be heated by microwave on low with 20 seconds. Prevented from violent boiling, the mixed reagents should be heated 10 seconds twice separately and waggled slightly in the heating gap.
2. Using the same protein with known concentration to prepare Standard Reagent can enhance the accuracy of concentration determining, e.g. determine the concentration of serum protein with Standard Reagent of BSA, or determine the concentration of antibodies with Standard Reagent of IgG.

Note:

1. Reagent A is clear liquid in the color of light yellow. Discard Reagent A that shows discoloration or sediments.
2. Reagents of BCA testing is slight toxic and alkaline. Please wear lab coat and gloves during assays.
3. BCA Protein Assay kit is intolerant with reducing or metal chelating reagents. If these reagents are inevitable, FineTest product- Bradford Protein Assay kit is available(**K002**)

Wuhan Fine Biotech Co., Ltd.

B9 Bld, High-Tech Medical Devices Park, No. 818 Gaoxin Ave. East Lake High-Tech Development Zone, Wuhan, Hubei, China(430206)

Tel: (0086)027-87384275

Fax: (0086)027-87800889

www.fn-test.com